
**Operating Instructions
Bedienungsanleitung
Mode d'emploi
Manual de Instrucciones
Istruzioni d'uso**

METTLER TOLEDO

**METTLER TOLEDO
BT-P42 Printer
Drucker BT-P42
Imprimante BT-P42
Impresora BT-P42
Stampante BT-P42**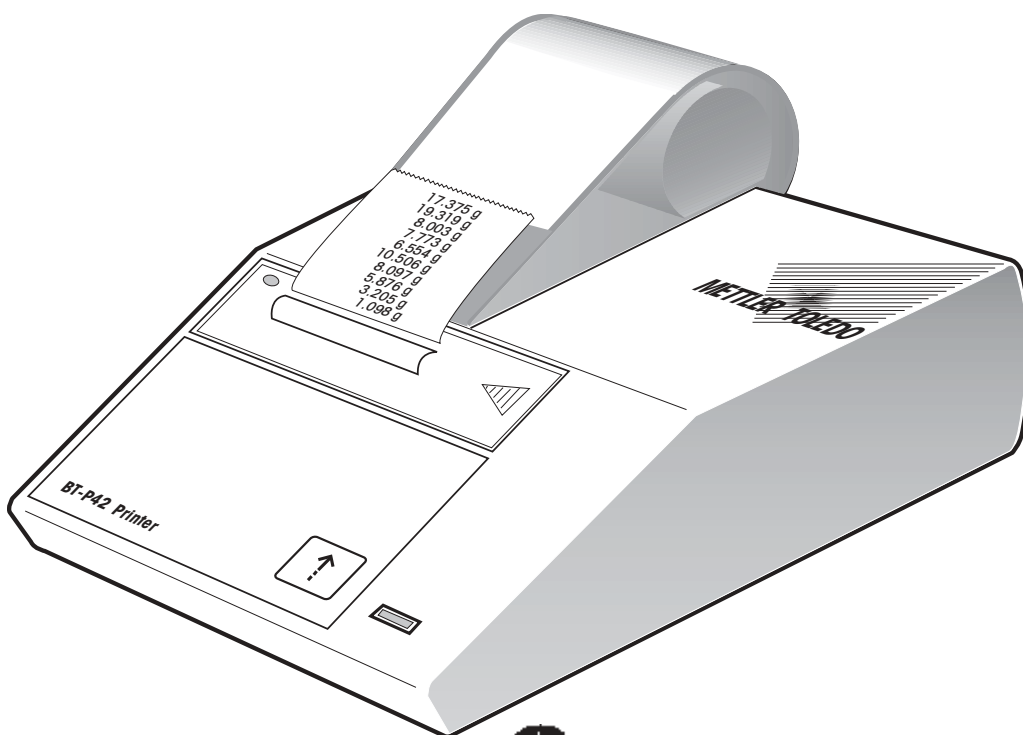

 **Bluetooth**® certified

Page 2

English

Seite 8

Deutsch

Page 14

Français

Página 20

Español

Pagina 26

Italiano

Content

1	Introduction	3
2	Startup procedure	3
2.1	Safety measures	3
2.2	Inserting paper	4
2.3	Connecting the printer to the device	4
3	Operation	4
4	Printer test	5
5	Maintenance	5
5.1	Changing paper	5
5.2	Replacing the ribbon	5
5.3	Cleaning	6
6	Malfunctions	6
7	Consumables	6
8	Technical data	7

1 Introduction

What are the capabilities of the BT-P42 Printer

The BT-P42 is a rugged and simple to operate dot matrix printer for METTLER TOLEDO balances with an a wireless Bluetooth® interface. The printout on normal paper fulfills the requirements of modern quality assurance systems (GLP, GMP, ISO 9001, etc.). These define, among other things, the principles and duties for the user of measuring instruments or inspection, measuring and test equipment. The BT-P42 complies with the principle of traceability of measurements by precisely documenting the origin of data, as well as calibration procedures. It fulfills the obligation to archive measurement results by its thermally stable and lightfast print. On connection to METTLER TOLEDO balances, the BT-P42 prints all data sent by the balance without restriction, that is not only weight values and calibration procedures, but also the results of all applications, built into the balance.

2 Startup procedure


2.1 Safety measures

The BT-P42 has been tested for the connections and intended purposes documented in these Operating Instructions. However, this does not absolve you from the responsibility of performing your own tests of the product supplied by us regarding its suitability for the methods and purposes you intend to use it for. You should therefore observe the following safety measures.


- The BT-P42 Printer may be used only indoors and on no account in hazardous areas.
- The electronics of the BT-P42Printer are protected against the ingress of splashing water and dust. However, as the printer housing is not tight the printer should not be used in the presence of liquids.
- Comply with cleaning instructions (see item 5.3), protect power plug against wet conditions.
- Never open the printer housing. Should you ever have problems with your printer, please contact your responsible METTLER TOLEDO dealer.

2.2 Inserting paper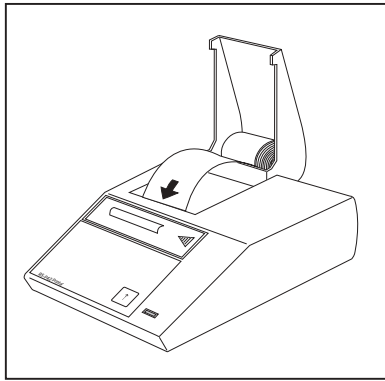

- Feed paper through the slot in the printing unit in the direction of the arrow.
- Switch on printer with toggle switch at the rear.
- Press and hold  until sufficient paper has been fed in.

2.3 Connecting the printer to the device

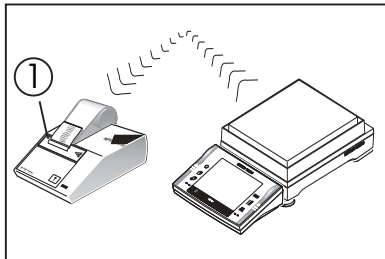- Position the printer within range of the balance.
Maximum distance: 10 m
- Connect the device in accordance with the Installation and Operating instructions for the BT or BTS interface option.

As soon as the device to be connected has been set up for the BT-P42, the printer is ready to operate. The connection may take 10 to 20 seconds to establish.

The connection status is indicated by the blue Bluetooth lamp (1):

- Lamp flashing: No connection
- Lamp continuously lit up: Connection present

3 Operation

- Pressing the printer key  advances the paper 1 line (line feed).
- Pressing the balance key with the  symbol prints out the stable weight value.
As long as the balance does not show a stable weight value, nothing will be printed out.

5.3 Cleaning

As the printer housing is made of top grade, resistant materials, all commercially available cleaning agents may be used.

6 Malfunctions

Malfunction / Message	Possible cause	Rectification
Printout unreadable	<ul style="list-style-type: none"> • Ribbon worn / entangled • Service life of printing unit at an end 	<ul style="list-style-type: none"> • Change / tension ribbon • Please contact METTLER TOLEDO service
Green pilot lamp does not light up	<ul style="list-style-type: none"> • No power • fuse defective 	<ul style="list-style-type: none"> • Printer on • Please contact METTLER TOLEDO service
Blue Bluetooth display does not light up	<ul style="list-style-type: none"> • No power • Bluetooth module faulty 	<ul style="list-style-type: none"> • Printer on • Please contact METTLER TOLEDO service
Blue Bluetooth display flashes continuously	<ul style="list-style-type: none"> • No connection • Bluetooth module faulty 	<ul style="list-style-type: none"> • Establish connection • Reduce distance • Please contact METTLER TOLEDO service

7 Consumables

	Type / Order number	Standard
Paper roll, set of 5	72456	2 piece
Paper roll, adhesive, set of 3	11600388	1 piece
Ribbon cartridge, black, minimum order quantity: 2	65975	1 piece
Spare printing unit (may be changed only by METTLER TOLEDO service!)	600133	—
Power line fuse 230 V (T100L250V) or 115 V (T200L250V) (may be changed only by METTLER TOLEDO service!)	commercially available	

8 Technical data

Printer

Printing unit	Dot matrix printer 5x9 dots, 24-character line length, Epson IBM character table No. 4
Printing speed	1 line per second
Ribbon cartridge	Exchangeable, black
Paper roll	Normal paper, 58 x dia. 51 mm, integrated in housing, commercial size

Printer functions

Printing, lightfast and thermally stable (GLP, GMP, ISO 9001), line feed, self-test

Data transmission

Wireless Bluetooth® technology
Max. range of 10 m

Designed test conditions

Power line voltage, frequency	115 V or 230 V, admissible fluctuations -20% +15%, 50/60 Hz
Power consumption	140 mA or 70 mA
Altitude	4000 m
Temperature range	0...40 °C
Humidity	Max. relative humidity 80% for temperatures to max. 31 °C, linear decreasing to 50% at 40 °C
Installation category	II
Pollution degree	2
Dimensions / Weight	W x D x H = 157 x 210 x 85 mm / net 1.5 kg (incl. paper roll)

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**To protect your METTLER TOLEDO product's future:
METTLER TOLEDO Service assures the quality, measuring accuracy and preservation of value of all METTLER TOLEDO products for years to come.
Please send for full details about our attractive terms of service.
Thank you.**

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P11780625

Subject to technical changes and to the availability of the accessories supplied with the instruments.