


Always on the ready the dual titration stand DV92

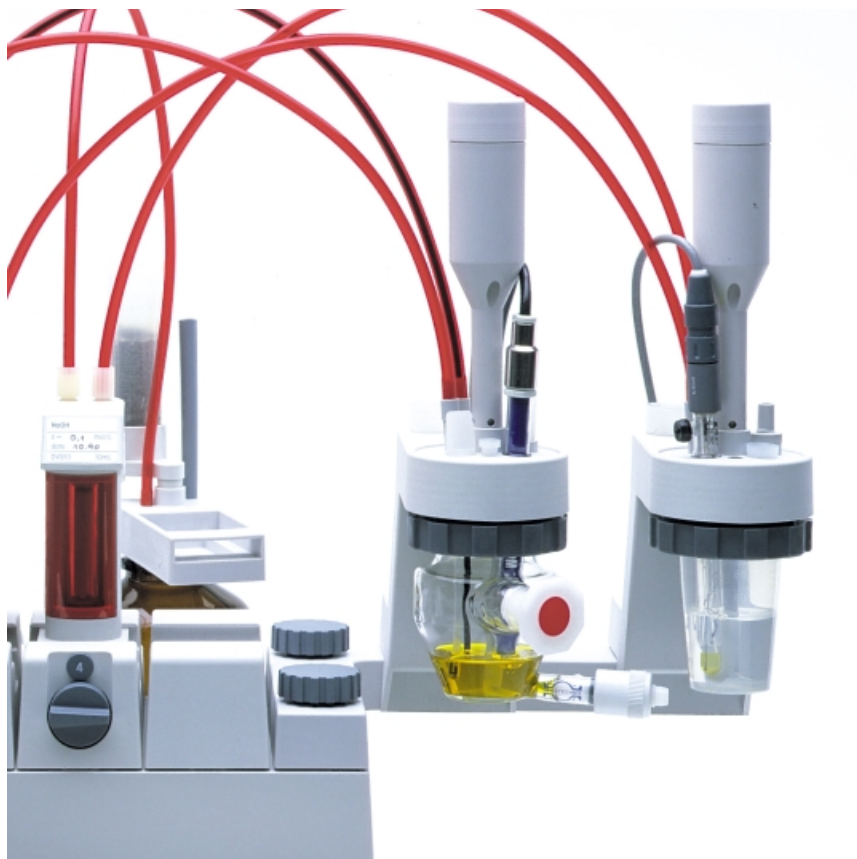

For many users it is a requirement that aqueous and non-aqueous titrations be performed on the same titrator. The changing over of sensors, burette tubes or stirrer can be precious time wasted in the laboratory.

By using the DV92 dual titration stand for the DL5x and DL7x family of titrators, it is possible to have two titration stands ready for use at all times. These stands can be arranged to accommodate aqueous titration on one and non-aqueous on the other. With the use of the optional DV750 KF set it is possible to have Karl Fischer titration of water content on the one stand and all other analyses on the other.

In many cases, where only two different analyses are performed on the titrator, it can simplify the analyst's life considerably by having one method set up per titration stand. In this way no changeover of titrants or sensors is required. The analyst simply attaches the sample to the appropriate titration stand and starts the method. The titrator knows from the method which stand is being used and automatically starts the correct stirrer. With the DL77 it is even possible to use both titration stands simultaneously thanks to its parallel titration capabilities.

The DV92 dual titration stand is delivered standard with an L-shaped arm to accommodate two titration stands, a second titration stand with accessories, and a second propeller stirrer.

Ordering information

Description	Part No.
DV92 dual titration stand	DV92
Optional for Karl Fischer titration:	
DV750 set for KF titration	23614
KF option for DL5x titrators	51107269
DK102A polarizing current source for DL7x titrators	106013
Double platinum pin electrode for KF with DK102A	DM142

subject to technical changes
© 12/99 Mettler-Toledo GmbH
Printed in Switzerland 51724110

Mettler-Toledo GmbH Analytical
Sonnenbergstrasse 74
CH-8603 Schwerzenbach
Tel. +41 1 806 77 11
Fax +41 1 806 73 50
<http://www.titration.net>

METTLER TOLEDO